



## Laboratory Guidebook Notice of Change

Chapter new, **revised**, or archived: MLG 5A.02

Title: FSIS Procedure for the Use of *Escherichia coli* O157:H7 Screening Tests

Effective Date: 10/01/10

Description and purpose of change(s):

Sections 5A.2 *Safety Precautions* and 5A.6.2 *Interpretation of Results* were clarified and Section 5A.1.2 *Limits of Detection* was removed. Raw beef mixed with other meat or poultry product was validated as an additional matrix for *E. coli* O157:H7 analysis.

**The methods described in this guidebook are for use by the FSIS laboratories. FSIS does not specifically endorse any of the mentioned test products and acknowledges that equivalent products may be available for laboratory use.**

**United States Department of Agriculture  
Food Safety And Inspection Service, Office of Public Health Science**

MLG 5A.02	Page 1 of 4	
Title: FSIS Procedure for the Use of <i>Escherichia coli</i> O157:H7 Screening Tests		
Revision: 02	Replaces: 01	Effective: 10/01/10

**Procedure Outline**

- 5A.1 Introduction
- 5A.2 Safety Precautions
- 5A.3 Quality Control Procedures
- 5A.4 Equipment, Reagents and Media
- 5A.5 Sample Preparation and Primary Enrichment
- 5A.6 The BAX<sup>®</sup> System and BAX<sup>®</sup> *E. coli* O157:H7 MP Test for Screening *Escherichia coli* O157:H7
  - 5A.6.1 Procedure
  - 5A.6.2 Interpretation of Results
- 5A.7 Use of Lateral Flow Devices for Screening *Escherichia coli* O157:H7
  - 5A.7.1 Test Procedure
  - 5A.7.2 Interpretation of Results
- 5A.8 Selected References

**5A.1 Introduction**

This method describes the use of a commercial PCR-based screening procedure to screen test enrichment cultures of ground beef, ground beef mixed with other meat or poultry product, fermented sausage, cooked meat patties, beef trim products, and environmental swabs for *Escherichia coli* O157:H7 (*E. coli* O157:H7). In the event that the BAX<sup>®</sup> system equipment or BAX<sup>®</sup> *E. coli* O157:H7 MP test kits are unavailable, this method describes the use of lateral flow devices that FSIS may use to screen for *E. coli* O157:H7.

Samples identified as potentially positive for the presence of *E. coli* O157:H7 by any of these tests are subject to cultural confirmation as described in MLG 5.

Unless otherwise stated all measurements cited in this method have a tolerance of  $\pm 2\%$ .

**5A.2 Safety Precautions**

*E. coli* O157:H7 is a human pathogen with a low infectious dose. (Ingestion of 100 cells can cause disease.) The use of gloves and eye protection is mandatory for all post enrichment viable culture work. Work surfaces must be disinfected prior to and immediately after use. Laboratory personnel must abide by CDC guidelines for manipulating Biosafety Class II pathogens. A Class II laminar flow biosafety cabinet is recommended for activities with potential for producing aerosols of pathogens. All available Material Safety Data Sheets (MSDS) should be obtained from the manufacturer for the media, chemicals, reagents and microorganisms used in the analysis. The personnel who will handle the materials should read all MSDS sheets.

**United States Department of Agriculture  
Food Safety And Inspection Service, Office of Public Health Science**

MLG 5A.02	Page 2 of 4	
Title: FSIS Procedure for the Use of <i>Escherichia coli</i> O157:H7 Screening Tests		
Revision: 02	Replaces: 01	Effective: 10/01/10

**5A.3 Quality Control Procedures**

- a. All media must be pre-warmed to 18-35°C prior to use.
- b. *E. coli* O157:H7 strain 465-97, or equivalent, shall be used as the positive control. See MLG 5, Section 5.3c for the positive culture control procedures.
- c. *E. coli* ATCC strain 25922, or equivalent, shall be used as the negative control for bead capture and screen tests.
- d. Prepare at least one “blank” (incubated but un-inoculated pre-enrichment/ enrichment broth) to provide a sterility control for the process.

**5A.4 Equipment, Reagents and Media**

In addition to equipment, reagents and media used in analysis of samples as described in MLG 5, the following materials will be needed.

- a. PCR tube holder (Qualicon or equivalent)
- b. Cell lysis tube cooling block (Qualicon or equivalent) held at  $5 \pm 3^\circ\text{C}$
- c. PCR cooling block (Qualicon or equivalent) held at  $< -10^\circ\text{C}$
- d. Heating block set at  $37 \pm 2^\circ\text{C}$
- e. Heating block set at  $95 \pm 3^\circ\text{C}$
- f. Repeating pipettor to deliver  $200 \pm 20 \mu\text{L}$ , and sterile tips
- g. Pipettor to deliver  $5 \pm 1 \mu\text{L}$ , and sterile disposable filtered tips
- h. Pipettor to deliver  $150 \pm 15 \mu\text{L}$ , and sterile disposable filtered tips
- i. Eight-channel pipettor to deliver  $50 \pm 5 \mu\text{L}$ , and sterile disposable tips
- j. 12 X 75 mm Falcon 352063, or equivalent, tubes
- k. Cell lysis tubes and caps, cell lysis tube rack and box (Genemate 8 strip tubes, ISC Bioexpress, T-3120-5 or equivalent)
- l. Pipettor or pipettes to deliver 5 mL
- m. BAX<sup>®</sup> System PCR Assay for Screening *E. coli* O157:H7 MP kit (Qualicon # 17720673) held at  $5 \pm 3^\circ\text{C}$
- n. Optional for Lateral Flow Device Testing:
  - i. Transia<sup>™</sup> Card *E. coli* O157:H7 (Diffchamb # ECO157), RapidChek<sup>®</sup> Pathogen Screening Test Kit (Strategic Diagnostics, Inc. # 7000160), or equivalent
  - ii. Polypropylene tubes and sterile pipets
  - iii. Heating block set at 97-100°C, boiling water bath, or autoclave (isotherm cycle)

**5A.5 Sample Preparation and Primary Enrichment**

Perform sample preparation and pre-enrichment as described in MLG 5, Section 5.5.

**United States Department of Agriculture  
Food Safety And Inspection Service, Office of Public Health Science**

MLG 5A.02	Page 3 of 4	
Title: FSIS Procedure for the Use of <i>Escherichia coli</i> O157:H7 Screening Tests		
Revision: 02	Replaces: 01	Effective: 10/01/10

**5A.6 The BAX<sup>®</sup> System and BAX<sup>®</sup> E. coli O157:H7 MP Test for Screening *Escherichia coli* O157:H7**

**5A.6.1 Procedure**

Following pre-enrichment incubation, perform the rapid screen using 5 ul of modified TSB broth with novobiocin plus casaminoacids (mTSB+n) pre-enriched samples. Follow the current BAX<sup>®</sup> System User's Guide for preparing reagents, performing the remainder of the PCR test, and reading the results. The equipment must be set up, operated, and all records documented according to laboratory work instructions.

**5A.6.2 Interpretation of Results**

- a. Samples that test BAX<sup>®</sup>-negative shall be reported as negative. Cultural analysis shall continue as per MLG 5, Sections 5.6 and 5.7 for sample pre-enrichments that test BAX<sup>®</sup> - positive, indeterminate, or have an invalid result. Or, the laboratory may review the cause and perform a correction. Based on the findings, the laboratory may analyze the indeterminate or invalid result samples by:
  - repeating the BAX<sup>®</sup> analysis from the rack loading step
  - preparing new BAX<sup>®</sup> tubes and repeating the analysis or
  - screen testing with a lateral flow device (LFD) described in Section 5A.7.
  
- b. In analytical runs where the positive control tests BAX<sup>®</sup>-negative, indeterminate, or has a signal-error result, the entire batch of samples is affected and a review of the cause and a correction shall be performed. Based on the findings the laboratory may analyze the samples by:
  - repeating the BAX<sup>®</sup> analysis from the rack loading step
  - preparing new BAX<sup>®</sup> tubes and repeating the analysis
  - screen testing with a lateral flow device (LFD) described in Section 5A.7
  - preparing fresh samples from reserve tissues or
  - analyzing all of the samples culturally.

**5A.7 Use of Lateral Flow Devices for Screening *Escherichia coli* O157:H7**

In the event that the BAX<sup>®</sup> system equipment fails or BAX<sup>®</sup> E. coli O157:H7 MP test kits are unavailable, the following lateral flow devices may be used until the BAX<sup>®</sup> system is available to screen for *Escherichia coli* O157:H7. The laboratory should follow the manufacturer's guidelines or other validated laboratory procedures. **RapidChek<sup>®</sup>** and **Transia<sup>™</sup>** lateral flow devices have been validated for use in FSIS laboratories.

**United States Department of Agriculture  
Food Safety And Inspection Service, Office of Public Health Science**

MLG 5A.02	Page 4 of 4	
Title: FSIS Procedure for the Use of <i>Escherichia coli</i> O157:H7 Screening Tests		
Revision: 02	Replaces: 01	Effective: 10/01/10

**5A.7.1 Test Procedure**

- a. Remove the appropriate number of test devices from storage.
- b. Transfer 1-3 ml of the enrichment broth (from each Stomacher bag) into a polypropylene tube with a sterile pipette. For use with a heating block, use ~ 1 ml. Heat the capped tubes at 97-100°C for the time specified by the manufacturer.
- c. Follow the current manufacturers' instructions for performing the test.

**5A.7.2 Interpretation of Results**

Follow the current manufacturer's instructions for reading the test device results.

**5A.8 Selected References**

Centers for Disease Control and Prevention and National Institutes of Health (CDC/NIH). 2007. BioSafety in Microbiological and Biomedical Laboratories, 5<sup>th</sup> ed. U.S. Government Printing Office, Washington, D.C. also found at the CDC internet site.

BAX<sup>®</sup> System PCR Assay for Screening *E. coli* O157:H7 MP User Guide Supplement and Automated Detection for Bacterial Screening User Guide, Dupont Qualicon.

RapidChek<sup>®</sup> *E. coli* O157: (including H7) Test User Guide, Strategic Diagnostics Inc.

Transia<sup>™</sup> Card *E. coli* O157 Manufacturer's Insert, Diffchamb